

Seals tightly, quickly and universally:
PARAFILM® M Sealing Film.




PARAFILM® M

F I R S T C L A S S · B R A N D

**The practical sealing film for
your lab.**

Whether it's a beaker or an Erlenmeyer flask, a weighing bottle or a graduated cylinder – the stretchable PARAFILM® M Sealing Film always fits. It protects the sample against contamination or evaporation and prevents spillage from upset vessels. PARAFILM® M – an indispensable accessory for your lab.

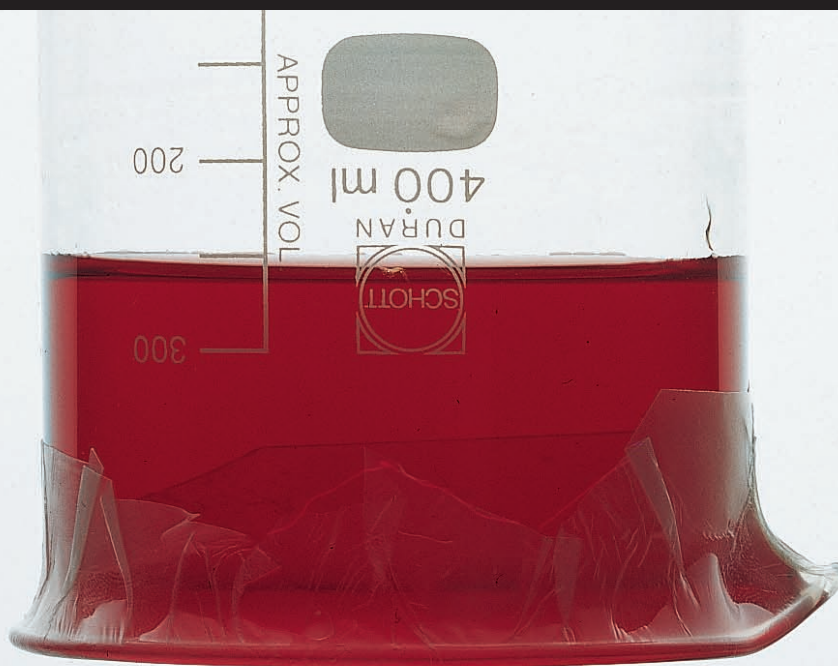

PARAFILM® M

Useful features

Simply cut PARAFILM® M off the roll and seal the vessel. After use, discard in an appropriate way.

PARAFILM® M always fits. It stretches up to 200% and clings even around irregular shapes and surfaces. Even a vessel falling over need not be a disaster. Provided it has been carefully sealed with PARAFILM® M, the contents will not spill out. (Vapor pressure from volatile substances may lift PARAFILM® M off the edge of the vessel.)



Physiological properties

PARAFILM® M is free of plasticizers, and consists largely of polyolefins and paraffin waxes. If PARAFILM® M should come into contact with food, the guidelines of the locally valid food handling laws should be followed.

Stretches to fit! PARAFILM® M Sealing Film

Chemical resistance

PARAFILM® M is resistant up to 48 hours against many polar substances, e.g., saline solutions, inorganic acids and alkaline solutions. After this period embrittlement may occur.

Storability

PARAFILM® M can be stored for at least 3 years without loss of quality under storage condition between 7 °C and 32 °C and a relative humidity of 50%.

Effects of 48 hours exposure at 23 °C:

Acids:	Hydrochloric acid 36.5%	resistant
	Sulphuric acid 98%	resistant
	Nitric acid 95%	resistant (brown coloration)
Alkaline solutions:	Sodium hydroxide 22%	resistant
	Ammonia 28%	resistant
Saline solutions:	Sodium chloride 20%	resistant
	Potass. permanganate 5%	resistant (brown coloration)
	Iodine solution 0.1 mol/l	resistant (brown coloration)
Organic solvents:	Methanol	resistant
	Ethanol	resistant
	Isopropanol	resistant
	Diethyl ether	not resistant
	Chloroform	not resistant
	Carbon tetrachloride	not resistant
	Benzene	not resistant
	Toluene	not resistant

PARAFILM® M complies with the general requirements of the FDA (Food and Drug Administration) for use under 55 °C, as well as satisfying GMP (Good Manufacturing Practice).

Cutter

PARAFILM® M Cutter

The handy PARAFILM® M Cutter is ideal for the neat and safe cutting of PARAFILM® M sealing film, and for its clean storage.



Very easy handling

The sturdy case of light-grey plastic stands safely on non-slip rubber pads. Two large turning wheels and smooth-running cutting blade make handling easy.

Comfortable and safe. PARAFILM® M Cutter



Inserting

Adjust the fitting bracket to the roll width (10 cm or 5 cm), and insert the PARAFILM® M carton with its flap folded over.



Dispensing

Turn the large wheels backwards until the desired length of PARAFILM® M appears.



Cutting

Cut off the adjusted length by moving the blade carriage across.



Taking

Another turn of the wheels releases the cut length.

Ordering Data

Technical Information

Physical properties

Toxicity: non-toxic
 Melting point: 60 °C
 Flash point: 301 °C
 Temperature range (continuous use): -45 °C to +50 °C
 Stretching ability: 200%
 Elongation at tear: 300%
 Gas permeability in 24 hours at 23 °C with 50% relative humidity:
 O_2 (oxygen): $\leq 350 \text{ cm}^3/\text{m}^2$
 N_2 (nitrogen): $\leq 105 \text{ cm}^3/\text{m}^2$
 CO_2 (carbon dioxide): $\leq 1100 \text{ cm}^3/\text{m}^2$
 Water vapor permeability in 24 hours at 37 °C and 90% relative humidity: $0.8 \text{ g}/\text{m}^2$



PARAFILM® M Cutter

Suitable for 50 mm and 100 mm wide rolls.
 Dimensions L x W x H:
 215 x 165 x 155 mm.
 Pack of 1.

Cat. No.	7016 50
----------	---------

Storability

PARAFILM® M can be stored for at least 3 years without loss of quality under storage condition between 7 °C and 32 °C and a relative humidity of 50%.

PARAFILM® M Sealing Film

Width mm	Length m	Pack of	Cat. No.
50	75	24	7016 11
100	38	12	7016 05
100	75	12	7016 06
500	15	6	7015 01

PARAFILM® M is available in four different sizes. The standard size is 100 mm wide and is supplied in rolls of 38 m or 75 m length.

BRAND® is a registered trademark of BRAND GMBH + CO KG, Germany.
 Other reproduced brands are the property of the respective owner.

Our technical literature is intended to inform and advise our customers. However, the validity of general empirical values, and of results obtained under test conditions, for specific applications depends on many factors beyond our control. Please appreciate, therefore, that no claims can be derived from our advice. The user is responsible for checking the appropriateness of the product for any particular application.

Subject to technical modification without notice. Errors excepted.

BRAND GMBH + CO KG · P.O. Box 11 55 · 97861 Wertheim · Germany
 Tel.: +49 9342 808-0 · Fax: +49 9342 808-236 · E-Mail: info@brand.de · Internet: www.brand.de

