

Pipetting comfort is just a PIVOT away



At a glance, monitor the conveniently top mounted speed control and low-battery indicator.

Rotate pipet nozzle to pipet in the most comfortable position possible.

Ergonomic design fits comfortably in your hand for all day use.

Concave pressure-sensitive trigger buttons provide precise control while reducing finger fatigue.

U.S. and European patents pending

Community Design No. 000203799-0001

Patent-pending adjustable nozzle feature gives you maximum pipetting control for a range of applications.



Position 1

Use in the standard vertical position for basic benchtop uses.



Position 2 — 7.5° Change

The smaller-angled position works well for tight benchtops with many bottles and flasks.



Position 3 — 15° Change

The wide-angle position is ideal for restricted spaces such as fume hoods or clean cabinets.

Rota-Filler 3000®

- Patent-pending adjustable nozzle rotates to three positions for maximum comfort and convenience.
- Contoured grip allows for relaxed hold to reduce hand strain.
- Three speed control options—high, low, and gravity


The Rota-Filler 3000® combines a powerful motor with a lightweight, comfortable design. Our commitment to the comfort of the end user lead us to make this an ergonomic filler. Features such as textured grip, hand-neutral design for right- and left-handed users, concave triggers, and convenient controls ensures the filler fits comfortably in your hand for many hours filling and dispensing liquid. The design has a patent pending as well.

Pipetting speed is easily set. Slide the speed switch between two operating modes—blow-out or gravity. Select the low or high range in "blow-out" mode to dispense the remaining liquid in the pipette. Flip to "gravity" mode for use with "To Contain" pipettes to fill in low mode and dispense the liquid without the motor.

Filler accommodates glass and plastic pipettes ranging in volume from 1 to 100 mL. The hydrophobic 0.45µm filter in the nozzle prevents contamination. The polypropylene pipet nozzle is autoclavable.

A red low-battery indicator signals when the batteries need to be recharged. Use filler during recharging, or place it in the included stand or holder. Filler includes two replacement 0.45µm filters, a dual-position charging stand, and a magnetic/wall-mountable holder. Filler measures 170 x 47 x 143mm.

Cat. No.	Description
HS3000RFAU	Rota-Filler 3000®, 110/120 VAC
HS3000RFAJ	Rota-Filler 3000®, 110/120 VAC
HS3000RFAT	Rota-Filler 3000®, 110/120 VAC
HS3000RFBC	Rota-Filler 3000®, 220/230 VAC
HS3000RFBA	Rota-Filler 3000®, 220/230 VAC
HS3000RFBI	Rota-Filler 3000®, 220/230 VAC
HS3000RFBK	Rota-Filler 3000®, 220/230 VAC
HS3000RFC	Replacement Filters, Pack of 5

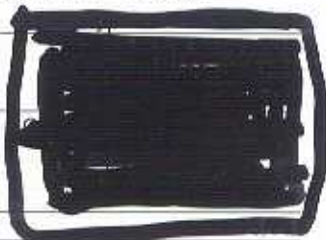

Dual-position stand conveniently charges the unit even with a pipette attached. Place unit upright or laying down.



Includes a handy magnetic wall mount holder—simply place on metal surface.

Holder can also be screwed in place on a non-magnetic surface.

Feedback From Customers



"I like the Rota-Filler 3000® better; I like being able to adjust the position of the pipet nozzle especially when working in confined spaces—like a clean cabinet or hood. The adjustable nozzle is a nice feature."

—M. Otegui, Assistant Professor

"The adjustable pipet nozzle is useful especially with tall, narrow bottles."

—A. Jacobs, Research Specialist



Manual Pipette Rack

Our FREE thank you gift for your purchase of the Rota-Filler 3000®. Holds 0.1 to 25 mL pipettes or cans. Two attachments hold Rota-Filler 3000® as well as pipet bulbs or pipetting aids. Measures 291 x 88 x 401mm.