



## Pipette filler, Bibbette®



- A novel PTFE and silicone rubber valve assembly controls the vacuum created by simply squeezing the unique rubber bulb
- By moving the thumb lever pipettes from 0.1ml up to 100ml can be easily and accurately filled and emptied
- Blow out pipettes are catered for by the depression of a small thumb operated squeeze ball
- A replaceable PTFE membrane filter ensures liquid is not accidentally drawn into the unit (2 supplied)
- The filter and pipette holding cone can be sterilised by autoclaving



Product Code	Description	Pack qty.
PFL001B	Bibbette pipette controller, Blue	1
PFL001N	Bibbette pipette controller, Plain	1

## Bibbette® accessories

Product Code	Description	Pack qty.
PFL002	Spare membrane filters pack of 10	10
PFL003	Spare suction bulb	1
PFL004	Spare valve system for Bibbette pipette controller	1
PFL006	Spare silicone adapter	1

## Pipette fillers, natural rubber



- Suitable for pipettes up to 1000ml
- Pipette filler works via a series of glass ball valves for air release, suction and delivery
- PWY145 is suitable for single handed operation
- PWY146 has an extended insert nozzle for improved grip

Product Code	Pack qty.
PWY144	1
PWY145	1
PWY146	1



### Operating instructions for PWY145

1. Evacuate via the automatic valve
2. Intake by pressure on operating point
3. Drain by pressure on operating point
4. Blow out the pipette (no valve actuation necessary)
5. Remove the valve for easy inside cleaning
6. Replace the valve using a twisting motion